Laser Diode Controllers

LDC500 series — Laser diode drivers with integrated temperature controllers



Laser Diode Driver

- · 100 mA, 500 mA or 2 A source
- Ultra-low drift (<10 ppm/°C)
- · <0.3 µA current noise (LDC500)
- · CC & CP mode dynamic switching

TEC Controller

- · 36 W output power
- Ultra-high stability (0.0005 °C/°C)
- · Thermistor, RTD and IC sensors
- · CC & CT mode dynamic switching



LDC500 Series Laser Diode Controllers

Introducing the LDC500 Series Laser Diode Controllers — highly stable, low-noise current sources, with integrated temperature controllers — all at very affordable prices.

The LDC500 series are the ideal instruments for controlling the current and temperature of your laser diodes. They have the performance and features you expect from instruments costing twice as much. The LDC501 has up to 500 mA of output current with less than 1.1 μA of rms noise, while the LDC500 has up to 100 mA of current and less than 0.3 μA of noise. The LDC502 provides up to 2 A of current.

With a low-noise current source, a 36 W high-precision temperature controller, and standard computer interfaces including Ethernet, the LDC500 series is the right choice for your laser diode testing and control applications.

Easy-To-Use Interface

The LDC500 series have an intuitive user interface, and many first time users will be able to operate the instrument without having to crack open the manual (although we do recommend reading the manual). Unlike competitive models, the LDC500 series controllers have a dedicated front-panel display for parameter entry. You don't have to sacrifice monitoring temperature or current to simply change an instrument setting — you have a separate two-line, blue alpha-numeric display for that. In addition, bright 5-digit green LED displays constantly monitor current and temperature, and are large enough to easily read from anywhere in the lab.



phone: (408)744-9040 www.thinkSRS.com

Laser Diode Protection

Multiple laser diode protection features, including slow start turn-on, adjustable current limits, and adjustable compliance voltage, keep your laser diodes safe when unexpected events occur. Fast clamping and shut down provide extra protection against intermittent contact with the laser. Combined, these features provide trouble-free, safe control of your laser diode.

Linear Power Supplies

Independent linear power supplies are used for the laser diode controller and the temperature controller. The supplies are designed with a magnetically shielded toroidal transformer, and provide ultra-clean, stable isolated power.

Computer Control

Remote operation of the LDC500 series is supported with GPIB, RS-232 and Ethernet interfaces. All instrument functions can be controlled and read over any of the interfaces. Up to nine complete instrument configurations can be saved in non-volatile RAM and recalled at any time. A TTL trigger output is also available to synchronize other test equipment.

Stable Laser Diode Controller

To ensure a stable optical output from your laser diode, the LDC500 series LD controllers were designed to deliver noise-free, precision operation. They are accurate to ± 0.01 % FS, have automated testing setups, noise is as low as 0.3 μA rms, and they meet a drift specification of 10 ppm/°C.

The controllers have two modes of operation for the laser diode: constant current and constant power. Constant current mode (CC) programs the source to a precise DC amplitude. Alternatively, the constant optical power mode (CP) servos the current source to maintain a constant signal on a monitor photodiode. Both control modes allow you to add an external

modulation signal, with adjustable bandwidth up to 1 MHz (in CC mode) or 10 kHz (in CP mode).

Another convenient feature that the LDC500 series offers is a fully programmable photodiode bias voltage. You can set the bias between 0 and 5 V from the front panel, or remotely using one of the computer interfaces.

Bumpless Transfer

A unique feature of the LDC500 series is dynamic "Bumpless Transfer" between CC and CP modes. This feature means you don't have to shut down your laser to switch modes — simply press the *Current/Power* button.

36 W Temperature Controller

The LDC500 series integrated 36 W temperature controller lets you adjust temperature with 0.001 °C resolution, and measure temperature with 0.01 °C accuracy (with a calibrated sensor). It maintains a typical stability of 0.0005 °C/°C with respect to room temperature, and has a very wide temperature control range.

The TEC controller also has two modes of operation: constant temperature mode (CT) controls the TEC current to maintain a fixed temperature (or raw sensor value), while constant current mode (CC) operates the TEC at a fixed current. Thermistor, RTD and IC sensors are all supported.

The LDC500 series offers an auto-tuning feature which automatically optimizes the PID loop parameters of the controller. Of course, full manual control is provided too. Dynamic transfer between CT and CC modes for the TEC is also easy — just press the *Temp/Current* button.



LDC500 front panel



LDC501 rear panel



Laser Diode Current Source

Current Source

Range 0 to 100 mA, 0 to 50 mA (LDC500)

0 to 500 mA, 0 to 250 mA (LDC501)

0 to 2 A, 0 to 1 A (LDC502)

Setpoint resolution 1 μA (LDC500), 10 μA (LDC501),

0.1 mA (LDC502)

Accuracy $\pm 0.02\%$ of full scale

Output impedance $>1 M\Omega$ (DC)

Stability

Thermal <10 ppm/°C Short-term (1 hr.) <5 ppm full scale Long-term (24 hr.) <15 ppm full scale

Noise (10 Hz to 1 MHz) LDC500

 $0.9\,\mu A \text{ rms}$ (high range / high BW) $0.6\,\mu A \text{ rms}$ (high range / low BW) $0.5\,\mu A \text{ rms}$ (low range / high BW) $0.3\,\mu A \text{ rms}$ (low range / low BW)

LDC501

4.5 μA rms (high range / high BW) 1.5 μA rms (high range / low BW) 2.3 μA rms (low range / high BW) 1.0 μA rms (low range / low BW)

LDC502

 $25 \,\mu A \, rms$ (high range / high BW) $5.0 \,\mu A \, rms$ (high range / low BW) $10 \,\mu A \, rms$ (low range / high BW) $3.5 \,\mu A \, rms$ (low range / low BW)

Compliance voltage

Range 0 to 10 V, programmable

Resolution 10 mV Accuracy 0.2 V

Current Limit

Resolution

Range 0 to 100 mA, 0 to 50 mA (LDC500)

0 to 500 mA, 0 to 250 mA (LDC501)

0 to 1 A, 0 to 2 A (LDC502) 10 μA (LDC500 & LDC501),

0.1 mA (LDC502)

Accuracy $\pm 0.1 \,\mathrm{mA}$ (LDC500 & LDC501),

±0.4 mA (LDC502)

Analog Modulation

Input range
Input impedance

 $-10\,\mathrm{V}$ to $+10\,\mathrm{V}$ $2\,\mathrm{k}\Omega$, typ.

Gain

CC mode LDC500

10 mA/V (high range) 5 mA/V (low range)

LDC501

50 mA/V (high range) 25 mA/V (low range)

LDC502

200 mA/V (high range) 100 mA/V (low range) CP mode LDC500 & LDC501: 500 μA/V

(PD current) LDC502: $1000 \,\mu\text{A/V}$

Bandwidth (-3 dB)

CC mode DC to 1.0 MHz

(high BW, LDC500 & LDC501)

DC to 0.8 MHz (high BW, LDC502)

DC to 10 kHz (low BW, all LDCs)

CP mode DC to 5 kHz (high bandwidth)

DC to 100 Hz (low bandwidth)

Monitor Photodiode

Bias voltage 0 to 5 V, programmable

PD current range 0 to 5,000 µA (LDC500 & LDC501)

0 to 10,000 μA (LDC502)

Setpoint resolution 0.1 µA (CP mode)

Setpoint accuracy $\pm 2 \,\mu\text{A} \,(\text{LDC500} \,\&\, \text{LDC501})$

 $\pm 4 \,\mu A \,(LDC502)$

Measurement & Display

Output current

Resolution $1 \mu A \text{ (LDC500)}, 10 \mu A \text{ (LDC501)},$

0.1 mA (LDC502)

Accuracy $\pm 0.02 \%$ FS

Photodiode current

 $\begin{array}{ll} Resolution & 0.1\,\mu A \\ Accuracy & \pm 0.02\,\%\;FS \end{array}$

Laser diode forward voltage

Resolution 1 mV

Accuracy $\pm 0.02 \%$ FS (4 wire)

Temperature Controller

Temperature Control

Control range

IC sensor -55 °C to +150 °C

Resistor sensor -150 °C to +250 °C $(10 \Omega \text{ to } 500 \text{ k}\Omega)$

Setpoint resolution

Temperature $0.001 \,^{\circ}\text{C}$ Resistance $0.1 \,\Omega$

Setpoint accuracy

 $\begin{array}{lll} \mbox{Temperature} & \pm 0.01\,^{\circ}\mbox{C (sensor dependent)} \\ \mbox{Resistance} & 0.1\,\% \mbox{ of sensor resistance} \\ \mbox{Stability (typ.)} & (\mbox{using a } 10\,\mbox{k NTC thermistor)} \\ \mbox{Thermal} & 0.0005\,^{\circ}\mbox{C/}^{\circ}\mbox{C (vs. ambient)} \\ \end{array}$

 $\begin{array}{ccc} \text{Short-term (1 hr.)} & \pm 0.001 \, ^{\circ}\text{C} \\ \text{Long-term (24 hr.)} & \pm 0.002 \, ^{\circ}\text{C} \\ \text{Control algorithm} & \text{PID} \end{array}$

Auto-tuning Step response

TEC Output

Source type Linear, bipolar current source

Current range $-4.5 \,\mathrm{A}$ to $+4.5 \,\mathrm{A}$



phone: (408)744-9040 www.thinkSRS.com $\begin{array}{lll} \text{Setpoint resolution} & 1 \text{ mA} \\ \text{Setpoint accuracy} & \pm 5 \text{ mA} \\ \text{Max. power} & 36 \text{ W} \\ \text{Compliance voltage} & > 8 \text{ VDC} \\ \end{array}$

Current noise <0.1 mA rms @ 1 A output

<0.2 mA rms @ 4 A output

Current limits

Range 0 to 4.5 AAccuracy $\pm 5 \text{ mA}$

Temperature Sensors

Thermistors $10 \text{ to } 500 \text{ k}\Omega (10 \,\mu\text{A}, 100 \,\mu\text{A},$

1000 μA excitation)

RTD Pt-100, Pt-1000 (1 mA excitation)

IC voltage sensors LM335 and equivalent IC current sensors AD590 and equivalent

Measurement & Display

Temperature

Resolution 0.001 °C

Thermistor

Resolution 0.1Ω

Accuracy $\pm 0.2\% + 0.05 \Omega$

TEC current

Resolution 1 mAAccuracy $\pm 5 \text{ mA}$

TEC voltage

Resolution 1 mV

Accuracy $\pm 5 \,\mathrm{mV} \; (4 \,\mathrm{wire})$

General

Instrument connectors DB9-F (laser diode), DB15-F (TEC)

BNC (modulate, trigger output)

Remote interfaces GPIB (IEEE488.2), RS-232, Ethernet

Power 100/120 VAC or 220/240 VAC,

 $50\,\text{Hz}/60\,\text{Hz},\,100\,\text{W}$

Dimensions $7" \times 5" \times 15"$ (WHL)

Weight 15 lbs.

Warranty One years parts and labor on defects

in material and workmanship

Ordering Information

LDC300	Laser Diode Controller
LDC501	Laser Diode Controller
LDC502	Laser Diode Controller
O500RM	Rack mount tray for LDC series
O500C1	6' LD cable w/ flying leads
O500C2	6' TEC cable w/ flying leads
O500C3	6' LD cable w/ D connector
O500C4	6' TEC cable w/ D connector
O500C5	12' LD cable w/ D connector
O500C6	12' TEC cable w/ D connector

About Thermal Stability

Temperature fluctuations in a typical laboratory environment can often exceed several degrees Celsius over the course of a day. Small temperature changes can mean significant current changes in your laser diode if your controller is not up to the task.

The LDC500, LDC501 and LDC502 have a temperature coefficient of 10 ppm/°C, which is a factor of five better than competing models, making it the ideal controller for precision laser diode experiments.

The graphs below demonstrate the temperature performance of the LDC500 series. Over a 24 hour period, the ambient temperature change in the laboratory exceeds two degrees Celsius. Note that the output current deviation of the LDC after warm-up is stable to better than ± 10 ppm.



